Digital scale CS3K

USER MANUAL

PRECAUTIONS

Never load the scale with more than the maximum capacity. Overloading will permanently damage it!

Always keep your scale in a clean environment. Dust, dirt, moisture, vibration, air currents and a close distance to other electronic equipment can all cause an adverse effect on the reliability and accuracy of your scale.

Avoid shaking, dropping, or subjecting the scale to shocks. This is a precision instrument and must be handled with extreme care.

Only operate the scale on a stable, vibration-free surface.

WEIGHING

1. Unlock for use by using the red switch located under the scale.
2. Place the scale on a horizontal flat surface, press [ON/OFF].
3. Wait until "0.0" is displayed.
4. Put the object(s) on the weighing platform.
5. Using the [UNIT] key, you can switch between the weighing units.

TARE WEIGHING

1. Turn on the scale as described above.
2. Place the tare item on the platform.
3. Press [ZERO] and wait until "0.0" is displayed.
4. Add the item to be weighed.

CALIBRATION

1. Turn on the scale by pressing [ON/OFF] key.
2. Wait until the LCD displays "0.0" and then press and hold [UNIT] key for 3 seconds. The LCD will display "CAL", and then release the [UNIT] key.
3. Press [UNIT] key again. The LCD will display "CAL" followed by the required calibration weight.
4. Add the required calibration weight. The display will show "PASS" and return to normal weighing mode.
5. Put away the calibration weight and press [ON/OFF] to turn off the scale and the calibration is finished.

TRANSPORTATION

During the transportation, please protect the loadcell by using the red switch located under the scale.

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